

Overview

- Uniquely engineered spherical microcarriers
- 150 – 300 μm in diameter
- Embedded paramagnetic particles manipulated by magnetic field
- Improved transfection
- Morphological control
- High survivability following cryopreservation
- Useful for high-throughput screening

Introduction

- Market pressure exists for new paradigms to meet the demands for more rapid and reproducible cell production to keep up with advances in cell-based assays used in drug discovery and development
- Novel automation platforms enable high productivity on three dimensional surfaces

Methods

- Global Eukaryotic Microcarrier (GEM)
- HeLa, HEK, and smooth muscle cells
- The Automated Partnership SelecT™
- New Brunswick Scientific Instruments NucleoCounter®
- Yellow Springs Instruments (YSI) 2700 SELECT™ Biochemistry Analyzer
- Nikon Eclipse T300 & TE 2000-E
- Kalypsys Systems high-throughput screening robot
- PerkinElmer® ViewLux™
- Amaxa Biosystems 96-Well Shuttle

An Emerging Revolution in Automated 3D Cell Culture

Cal Trepagnier (1,2), Stacey L. Szymanski (1), Kevin W. Huff (1), Alison Rush (1), Richard Peltier (1), Berta Strulovici (1), Rebecca C. Lobo (3), Robin A. Felder (2)

(1) Department of Automated Biotechnology, Merck Research Laboratories, North Wales, Pennsylvania, United States of America

(2) Medical Automation Research Center, University of Virginia, Charlottesville, Virginia, United States of America

(3) BioCell Cubed, Inc. (formerly Global Cell Solutions, LLC), Charlottesville, Virginia, United States of America



Acknowledgements: Raul Lacson (1), Eric Johnson (1), Jason Cassaday (1), Shi Xiaoqing (1), Tarak Shah (1)

Results

Transfection

- GFP expression was uniform and in similar concentrations in cells on GEMs as compared to conventional transfection (Figure 1)
- HeLa cells stayed on the microcarriers after the electroporation procedure

Morphological control

- 3D surface on GEMs resulted in even confluent growth (Figure 2) as compared to clustering effect on 2D plate
- Smooth muscle cells wrapped around GEMs (Figure 3A)
- HEK cells more columnar than those grown on a tissue culture flask (Figure 3B)

Cryopreservation

- HeLa cells were cryogenically preserved on microcarriers for three weeks
- Forskolin stimulated cAMP dose response showed similar functionality after thaw and recovery (Figure 4)

High-throughput screening

- GEMs covered with cells were readily dispensed from a Kalypsys Systems kinetic straight-tip bottle-fed dispensing robot into the wells (Figure 5)



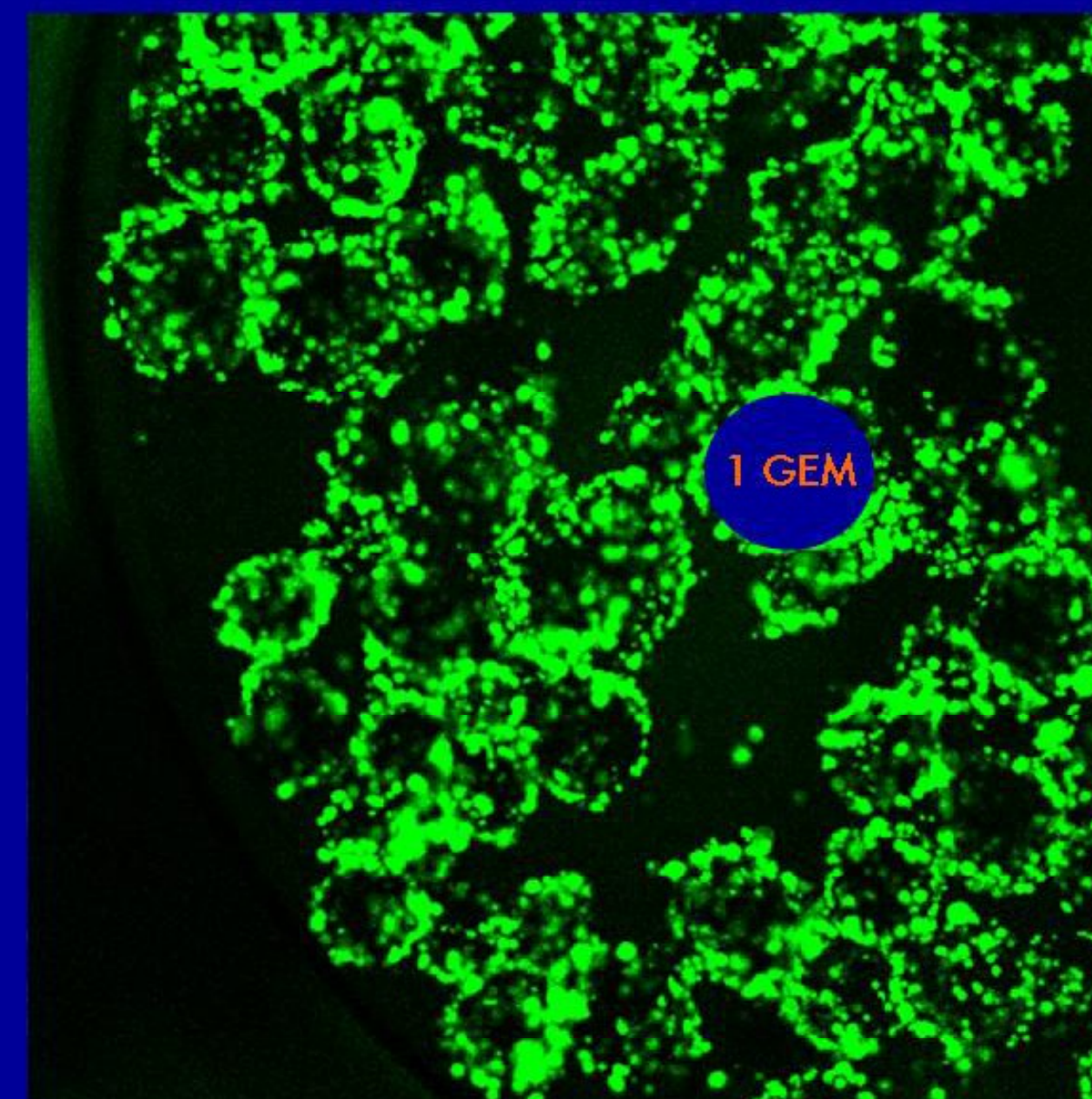


Figure 1 (above): Transfection of HeLa cells to express green fluorescent protein (GFP)
• Image taken at 24 hours using a confocal Nikon Edipse T300 with argon ion laser hookup at 40X and 7.40 gain

HeLa cells (100X)

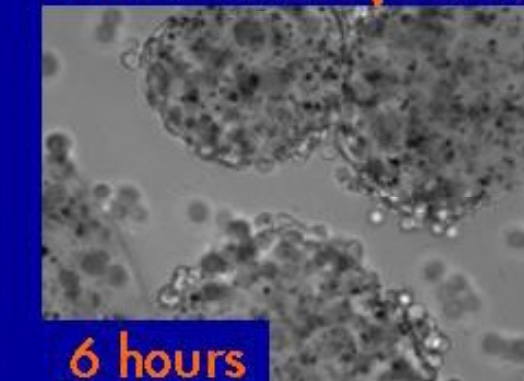


6 hours

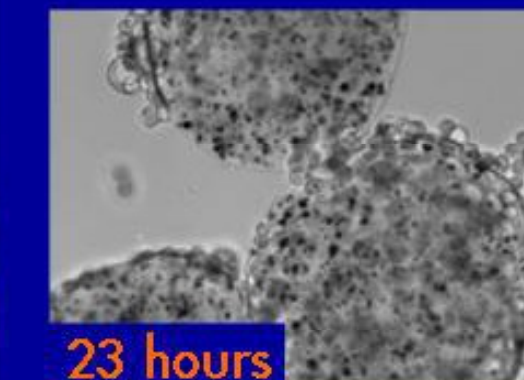

23 hours



48 hours

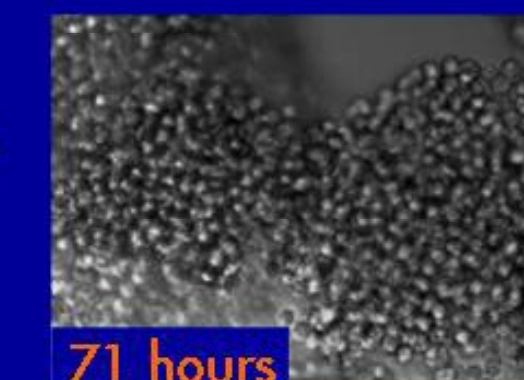

71 hours

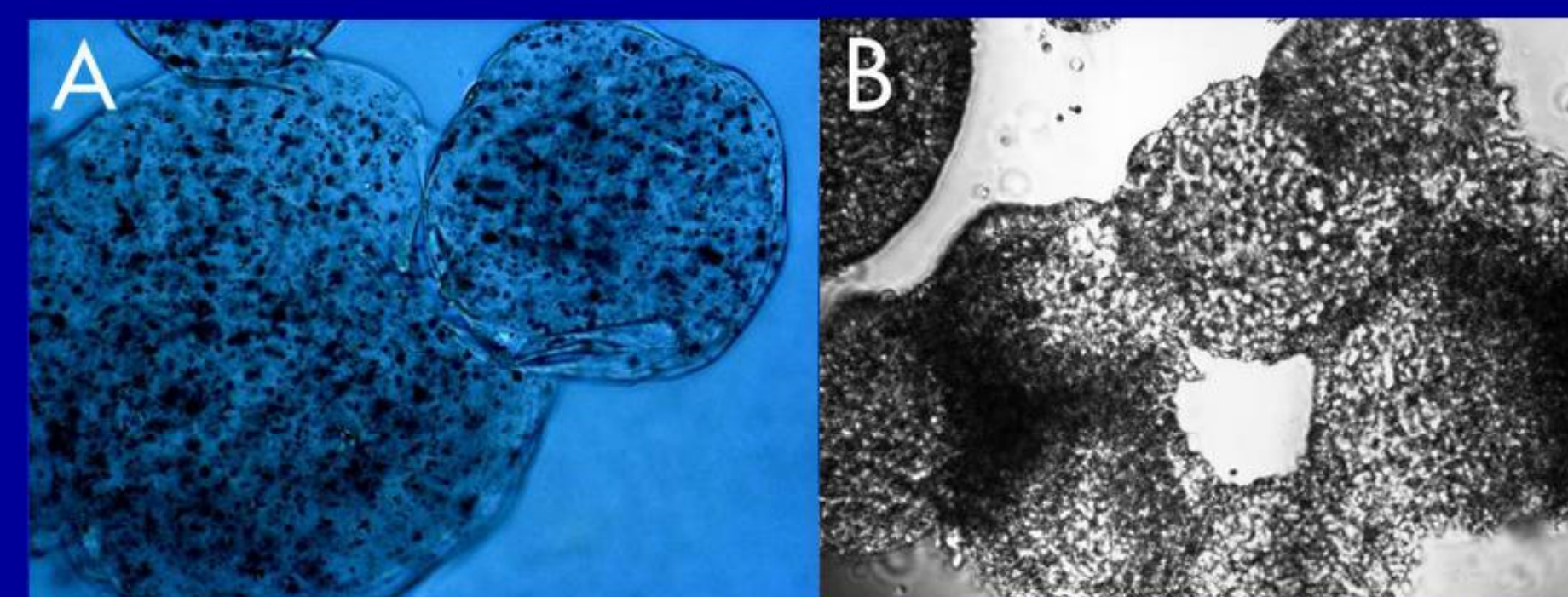


Figure 3 (above): Morphological control of microcarriers (A) smooth muscle cells at 24 hrs with blue filter (B) HEK cells at 5 days. Images taken with Nikon Edipse TE2000-E at 40X



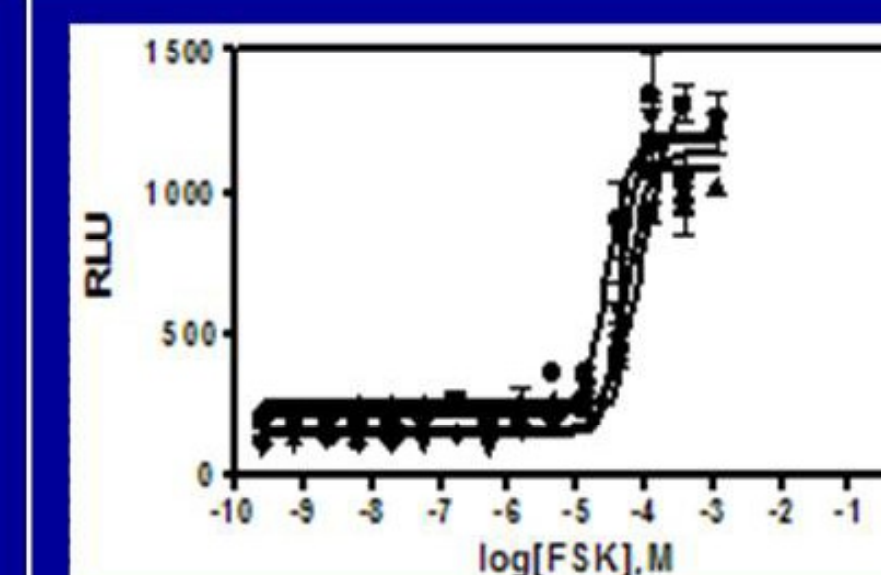


Figure 4:
Recovery following cryopreservation and pharmaceutical response
■ Cells frozen by traditional means
▲ Fresh cells
▼ Cells frozen on microcarriers at 50% confluence
● Cells frozen on microcarriers at 60% confluence
◆ Cells frozen on microcarriers at 90% confluence
No significant difference was observed between the five conditions

Figure 5:
High-throughput screening robotic dispense on to a 1536-well microplate
• The Cell Titer Glo® confirmation showed fluorescence that was measured by a PerkinElmer® ViewLux™
• Relative number of viable cells undergoing metabolism revealed by test for ATP

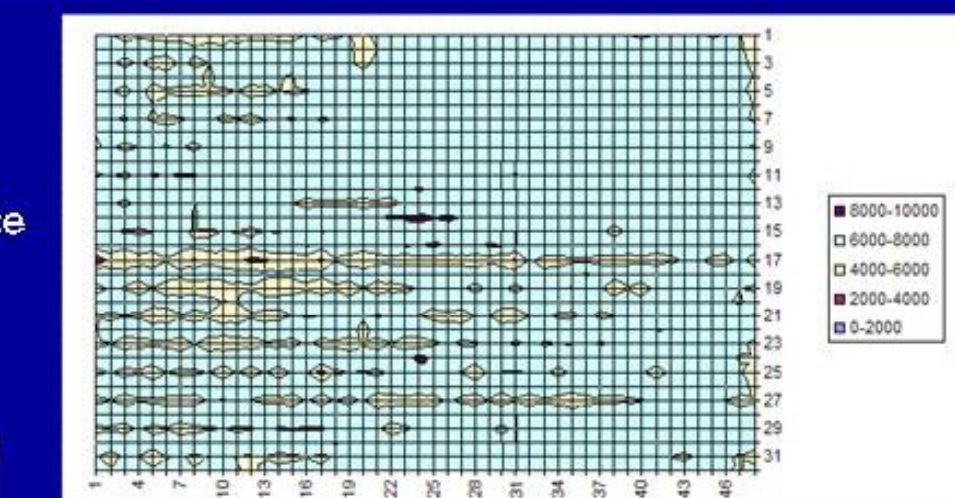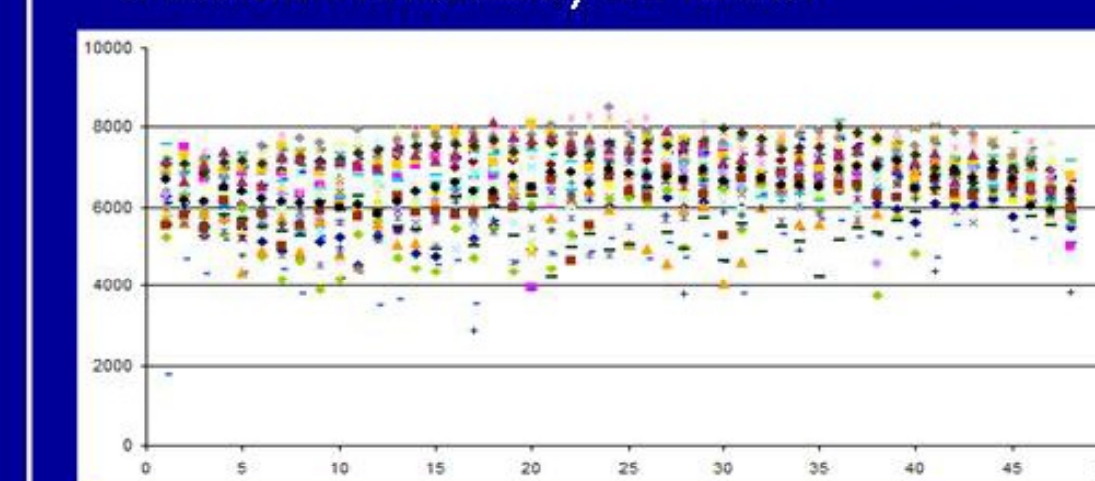

Vital statistics from heat map (above) and scatter plot (left):
• Mean: 6612.8 RLU/well
• Standard Deviation: 869.6 RLU/well
• Coefficient of Variation: 13.15 RLU/well
• Data suggests even cell distribution in plate

Conclusions

- Magnetic microcarrier culture offered improvements in yield, functionality, and convenience for all cells tested
- Cells grew to confluence in three days or fewer, averaging 750 cells/GEM
- Microcarrier-based cell manipulation was efficient for electroporation, robotic dispensing, and cryopreservation
- The convenience of microcarrier-based culture provided functional convenience of being able to pipette cells directly into a 1536-well microplate used in high-throughput screens without the need to passage